


Energy Transfer - Wind Turbine

Model No. ET-8783

Equipment List

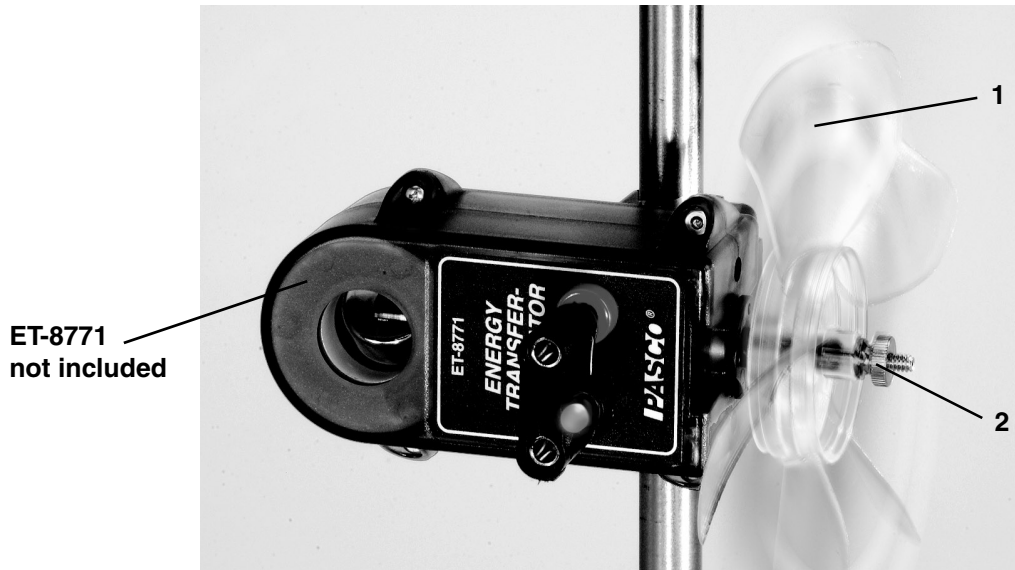

| Included Equipment | Replacement Model Number* |
|------------------------------|---------------------------|
| 1. 5-Blade Fan | 550-023 |
| 2. Thumb Nut | 614-006 |
| 3. Spacer (hidden from view) | 615-046 |

*Use Replacement Model Numbers to expedite replacement orders.

| Additional Equipment Required | |
|-------------------------------|---------|
| Energy Transfer Generator | ET-8771 |

Introduction

The Energy Transfer - Wind Turbine accessory can be attached to the Energy Transfer Generator (ET-8771) to create a wind energy turbine.

Equipment Setup

Energy Transfer-Wind Turbine Assembly

1. Remove the thumb nut and pulley from the Energy Transfer Generator. Do not discard these components. Retain for other use.
2. Place spacer over plastic shaft of the Energy Transfer-Generator.
3. Place 5-blade fan over same plastic shaft of Energy Transfer Generator.
4. Note orientation of 5-blade fan. See Figure 1.
5. Secure 5-blade fan with thumbnut. Do not over tighten.

