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### INFORMATION GENERALE

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### GENERAL INFORMATION

- 1) Handle the parcel with care. Unpack and check that the contents coincide with the packing-list. If any part is damaged or missing, please advise the distributor immediately.
- 2) Do not install or use the equipment without reading this handbook before.
- 3) This handbook must always be attached to the equipment and it must be available for all users.
- 4) If you have any doubts or enquiries, please contact with your supplier or J.P. Selecta.
- 5) **IMPORTANT! J.P. SELECTA WILL NOT ACCEPT ANY APPARATUS TO BE REPAIRED IF IT IS NOT DULY CLEANED.**
- 6) If any modification, elimination or lacking in maintenance of any device of the equipment by the user transgress the directive 89/655/CEE, the manufacturer is not responsible for the damage that can occur.
- 7) Do not use the apparatus with liquids which can give off vapours capable of making explosive mixtures.

### CONTENTS

|                             |
|-----------------------------|
| GENERAL INFORMATION         |
| CONTENTS                    |
| PACKING LISTS               |
| TECHNICAL FEATURES          |
| EQUIPMENT DESCRIPTION       |
| INSTALLATION                |
| OPERATION                   |
| MAINTENANCE                 |
| ELECTRICAL DIAGRAM          |
| GUARANTEE                   |
| "EC" CONFORMITY DECLARATION |

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### PACKING LIST

The standard equipment consist of the following components:

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## TECHNICAL FEATURES

Voltage supply 115-230V 50/60 Hz according to the characteristics plate indications.

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## EQUIPMENT DESCRIPTION

CENCOM is a benchtop centrifuge with an angle head of 6 15ml. tubes. It is supplied with safety cover and 8 speed positions.

### INSTALLATION

Place the centrifuge on a flat, horizontal, level surface.  
Connect the cable into the mains.

**ATTENTION! IMPORTANT POUR VOTRE SECURITE  
CAUTION!!! IMPORTANT FOR YOUR SAFETY**

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Do not use the apparatus to centrifuge liquids which are susceptible to form explosive mixtures.

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Be sure that the voltage supply is the same as the one indicated on the characteristics plate of the apparatus.

Do not use the apparatus if it is NOT EARTHED.

If you have to change the plug bear in mind the following:

Blue cable: Neutral.  
Brown cable: Phase.  
Yellow/green cable: Earth.

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### STARTING UP:

1. Connect into mains.
2. Put the tubes into the head in a balance way. If all buckets are not used, take notice that the load must be balanced symmetrically.
3. Close the cover.
4. Select the speed. (r.p.m.): 1000, 1500, 2000, 2500, 3000, 3500, 4000.
5. Select the centrifugation time. (hh.mm)
6. Press On / Off to start centrifugation.

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### STOP:

Press **On / Off** and wait few minutes before opening the door.  
 Do not place tubes and remove if the head is not complete stopped.  
 Leave the head to stop by itself. Do not use any tool or your hand to brake the head.

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### SAFETY:

Microswitch which disconnects the motor when opening the cover (in case of stop omission of the motor)

**CAUTION!! DO NOT OPEN THE COVER WHEN THE MOTOR IS WORKING**

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### MAINTENANCE

Before removing the cover disconnect the apparatus from the mains.

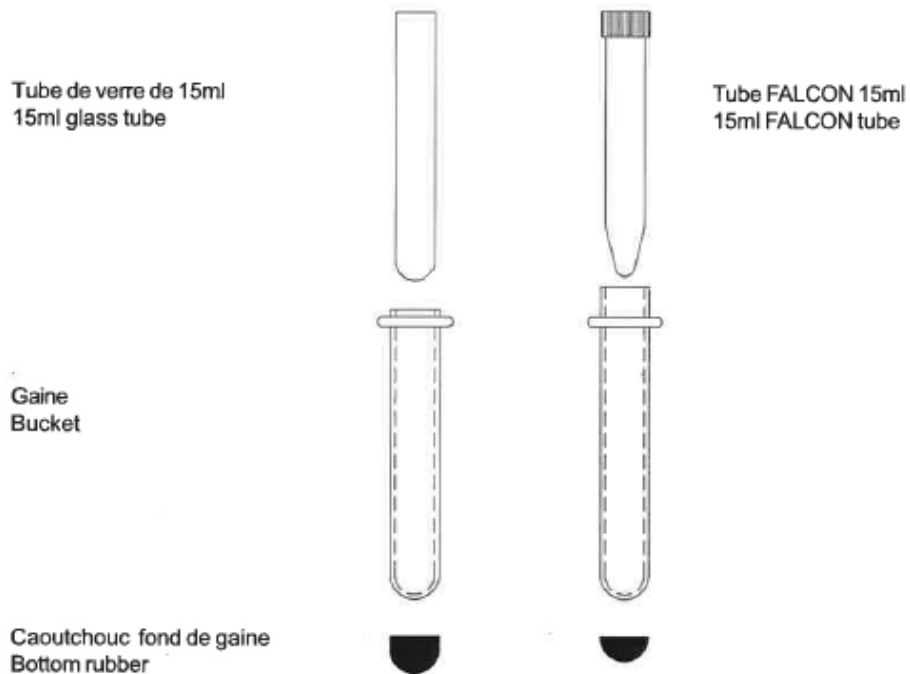
#### CLEANING:

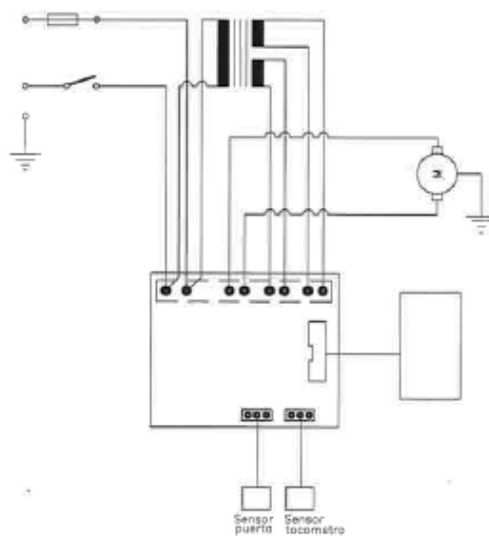
For the cleaning of the different parts use the following products:  
 Cleaning of casing, plastic and painted metals: Alcohol with cotton duster.

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